



## Installing a new IMGT library in SBTengine

1. Download the IMGT from the GenDx website (you can find the libraries following this link: <https://www.gendx.com/database-downloads-SBTengine/>)
2. Save the IMGT at this location: "C:\ProgramData\GenDx\Shared\Libs\UserLibs"
3. Start SBTengine and go to Help -> Preferences. Click on 'Library and Data folders'.
4. At 'Allele libraries', select Custom, then select the latest IMGT from the dropdown:

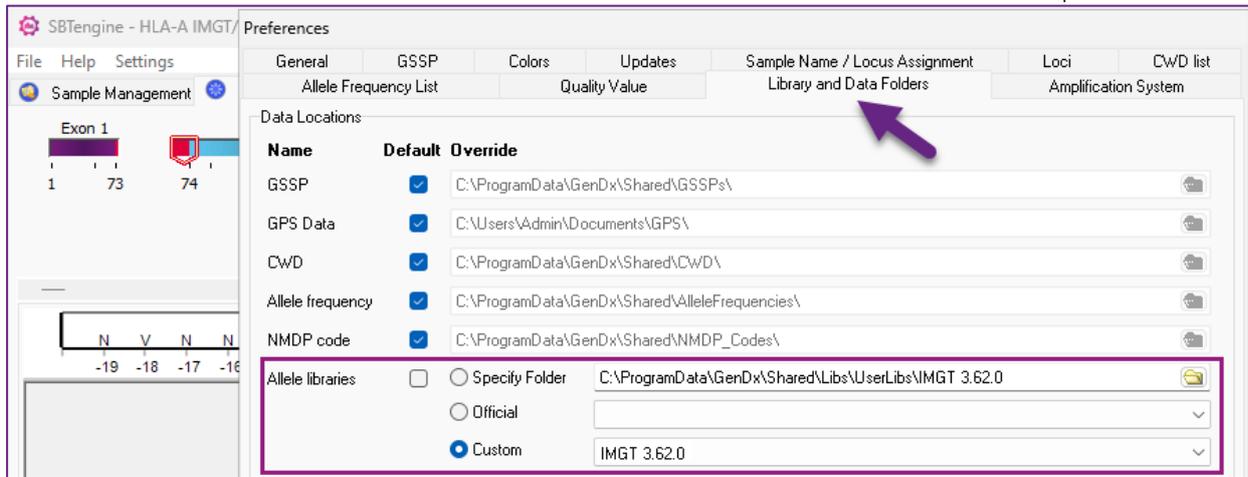

5. Select OK to save. After a few seconds, the new IMGT is shown in the header of the software:

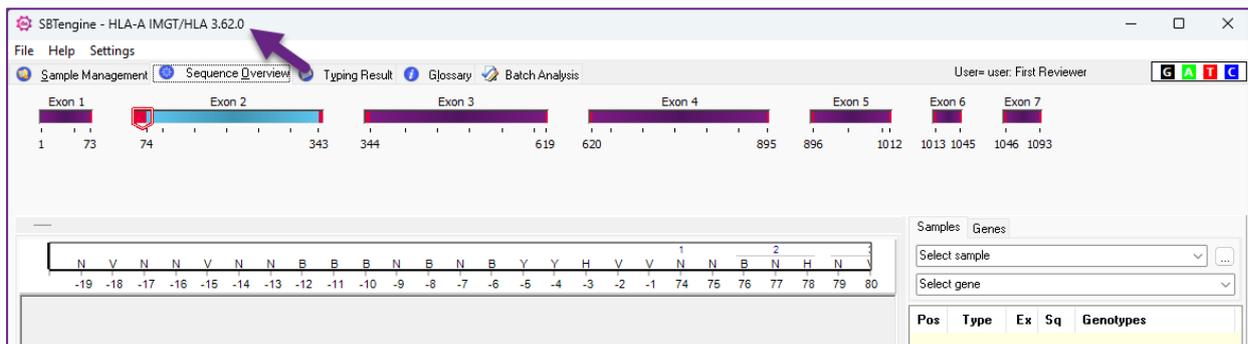

6. SBTengine now supports secure analysis of Sanger data with the latest IMGT/HLA Database.